

UiO : Faculty of Medicine		
Guidelines for use of the Saveen transilluminators and the BiO RAD Doc XRS geldokumentation		Version 2: 26.05.16 Version : 1 - 2009
		Page : 1
Elaborated for NCMM : HSE coordinator: Liv Bjørland	Approved:	Date : 25.05.16, Converted: 17.02.17

Saveen transilluminator: T-15M 312nm

BIO RAD transilluminator with trans-UVB 302nm

The ultraviolet transilluminators are a powerful source of UV-radiation that will cause damage to unprotected eyes and skin. Before operating the units be sure you are protected if the UVB-source is used.

When exposed:

- A UV- face mask must to be used when the light cover plate is removed / not in use.
- Gloves and long sleeves must be used to cover arms and hands to avoid sun burns.
- If another UV- source is installed please inform the people in charge and the safety delegate.