

Instructions for UVC-disinfection.

Version 2: 09.12.08

Version : 3

Page : 1 of 1

Elaborated for NCMM :

Earlier Radiation safety coordinator: Jorun

Solheim. Safety deputy: Gladys Tjørhom

HSE coordinator: Liv Bjørland

Approved:

Kjetil Taskén

Date : 17.02.15

Converted:17.02.17

Instructions for use of UVC in the ceiling in room 52:059, E3

- It is not allowed to stay in the room when the UVC-light in the ceiling is activated.
- The use of UVC for disinfection of the room is no longer routine, and was previously always arranged with the radiation safety coordinator.
- During the use of UVC the room must be labeled: **DANGER UVC**.

To avoid any accidental use of the UVC-light by the cleaning personnel, the fuse has been inactivated.

Instructions for use of UVC in the LAF-benches.

- The use of UVC- disinfection is not recommended, but can be used briefly either before or after work in the bench.
- Never leave the UVC on for a long time. The UVC will affect the equipment in the bench.
- The hinged window in front will protect the person from the UVC, if in use
- The UVC will not switch on if the hinged window is not properly closed.